

Detection of WNT2 in Formalin-Fixed, Paraffin-Embedded Mouse Tissue

Reagent and Antibody Information

[1X Wash Buffer](#)

[3% Hydrogen Peroxide](#)

[1% BSA Diluent](#)

[1X Citrate Buffer](#)

[Normal Rabbit IgG – Affinity Purified](#)

[DAB Chromogen](#)

[Hematoxylin](#)

Blocking Serum: Normal Donkey Serum

Jackson ImmunoResearch Laboratories, Inc.

West Grove, PA 19390

www.jacksonimmuno.com

1-800-367-5296

Catalog # 017-000-011

Avidin / Biotin Blocking Kit

Vector Laboratories, Inc.

Burlingame, CA 94010

www.vectorlabs.com

1-800-227-6666

Catalog # SP-2001

Primary Antibody: Rabbit Monoclonal WNT2 Antibody

Epitomics, Inc.

Burlingame, CA 94010

www.epitomics.com

1-877-772-2622

Catalog # 3169-1

Secondary Antibody: Donkey Anti-Rabbit IgG (H+L) Biotin-SP-Conjugated

Jackson ImmunoResearch Laboratories, Inc.

West Grove, PA 19390

www.jacksonimmuno.com

1-800-367-5296

Catalog # 711-065-152

Label Complex: R.T.U. Vectastain Elite ABC Reagent

Vector Laboratories, Inc.

Burlingame, CA 94010

www.vectorlabs.com

1-800-227-6666

Catalog # PK-7100

Staining Procedure

Positive Control Tissue: Female reproductive tract or papilloma skin

Stain Localization: Cytoplasmic and membrane

1. Deparaffinize and hydrate slides through the following solutions:

Solution	Repetitions	Time
Xylene	2 times	5 minutes
100% Ethanol	2 times	3 minutes
95% Ethanol	2 times	3 minutes
1X Wash Buffer	2 times	5 minutes

2. Quench endogenous peroxidase by placing the slides in 3% hydrogen peroxide for 15 minutes.

3. Rinse slides in 2 changes of 1X wash buffer for 5 minutes each.

4. Heat-Induced Epitope Retrieval Using The Steamer

Add distilled water to the bottom portion of the steamer.

Preheat 200ml of 1X citrate buffer in a Tissue Tek® container in the steamer between 95°C and 100°C.

Immerse a full rack of slides into the citrate buffer and place the lid back on the steamer.

(Insert blank slides into any empty slots in the rack to ensure even heating of slides)

Steam the slides for 20 minutes.

Remove container from steamer and cool for 20 minutes.

Rinse the slides in 2 changes of distilled water for 3 minutes each time.

5. Rinse slides in 2 changes of 1X wash buffer for 5 minutes each.

6. Block with 10% normal donkey serum for 20 minutes at room temperature.

Lot # _____ Date Reconstituted _____

DO NOT RINSE SLIDES. CONTINUE TO AVIDIN-BIOTIN BLOCK.

7. Avidin / Biotin Blocking Kit

Lot # _____ Exp. Date _____ New Kit: yes / no

Apply avidin block for 15 minutes at room temperature.

Quick rinse in 1X wash buffer.

Apply biotin block for 15 minutes at room temperature.

DO NOT RINSE SECTIONS WITH BUFFER BEFORE ADDING PRIMARY ANTIBODY.

ONLY WIPE EXCESS BLOCK.

8. Apply primary antibody at a 1:500 dilution. Incubate for 1 hour at room temperature.

Lot # _____ Exp. Date _____

For negative control slides, dilute normal rabbit IgG so that it's IgG protein concentration matches that of the primary antibody (if necessary). Then make a 1:500 dilution. If the concentrations can't be matched using this method, the dilution for the negative reagent may need to be adjusted. Apply the

negative and incubate for 1 hour at room temperature.

Lot # _____ Exp. Date _____

9. Rinse the slides in 2 changes of 1X wash buffer for 5 minutes each.

10. Apply the donkey anti-rabbit secondary antibody at a 1:1000 dilution. Incubate for 30 minutes at room temperature.

Lot # _____ Date Reconstituted _____

11. Rinse the slides in 2 changes of 1X wash buffer for 5 minutes each.

12. Apply the Vectastain R.T.U. Elite Label and incubate for 30 minutes at room temperature.

Exp. Date _____ New Kit: yes / no

13. Rinse the slides in 2 changes of 1X wash buffer for 5 minutes each time.

14. Apply the DAB chromogen. Incubate in the dark for 6 minutes at room temperature.

(Add 1 drop of DAB per ml of substrate)

Lot # _____ Exp. Date _____ New Kit: yes / no

15. Rinse the slides in tap water 3 minutes.

16. Counterstain with hematoxylin for 20 seconds.

17. Rinse the slides in tap water until water is clear.

18. Gently agitate slides in 1X wash buffer until the tissues turn blue.

19. Dehydrate through the following solutions:

Solutions	Repetitions	Time
95% Ethanol	1 time	3 minutes
100% Ethanol	3 times	3 minutes
Xylene	2 times	5 minutes

20. Coverslip

Updated 06/01/11