

Detection of Von Willebrand Factor (Factor 8) in Formalin-Fixed, Paraffin-Embedded Human Tissue

Reagent and Antibody Information

[1X Wash Buffer](#)
[3% Hydrogen Peroxide](#)
[1% BSA Diluent](#)
[Carezyme II \(Pepsin\)](#)
[DAB Chromogen](#)
[Hematoxylin](#)

Staining Kit: Vectastain Elite ABC Kit (Rabbit IgG)

Vector Laboratories, Inc.
Burlingame, CA 94010
www.vectorlabs.com
1-800-227-6666
Catalog # PK-6101

Note: This kit contains all reagents necessary to make the blocking reagent, secondary antibody and label complex.

Avidin / Biotin Blocking Kit

Vector Laboratories, Inc.
Burlingame, CA 94010
www.vectorlabs.com
1-800-227-6666
Catalog # SP-2001

Primary Antibody: Rabbit Anti-Human Von Willebrand Factor (Factor 8) Antibody

Biocare Medical
Concord, CA 94520
www.biocare.net
1-800-799-9499
Catalog # CP 039

Negative Control Serum: Normal Rabbit Serum

Jackson ImmunoResearch Laboratories, Inc.
West Grove, PA 19390
www.jacksonimmuno.com
1-800-367-5296
Catalog # 011-000-001

Staining Procedure

Positive Control Tissue: Endothelial cells
Stain Localization: Cytoplasmic

1. Deparaffinize and hydrate slides through the following solutions:

Solution	Repetitions	Time
Xylene	2 times	5 minutes
100% Ethanol	2 times	3 minutes
95% Ethanol	2 times	3 minutes
1X Wash Buffer	2 times	5 minutes

2. Quench endogenous peroxidase by placing the slides in 3% hydrogen peroxide for 15 minutes.
3. Rinse the slides in 2 changes of 1X wash buffer for 5 minutes each.
4. Proteolytic-Induced Epitope Retrieval Using Pepsin
Incubate the slides in Carezyme II: Pepsin (predilute) for 5 minutes at 37°C.
(Allow the pepsin to reach room temperature before use.)
Rinse the slide in distilled water for 1 minute to stop the enzymatic reaction.
5. Rinse the slides in 2 changes of 1X wash buffer for 5 minutes each.

Vectastain Rabbit Elite Staining Kit

Exp. Date _____ New Kit: yes / no

6. Apply the block from the Rabbit Elite Kit. Incubate for 20 minutes at room temperature.

DO NOT RINSE THE SLIDES. CONTINUE TO AVIDIN-BIOTIN BLOCK.

7. Avidin / Biotin Blocking Kit

Lot # _____ Exp. Date _____ New Kit: yes / no

Apply avidin block for 15 minutes at room temperature.

Quick rinse in 1X wash buffer.

Apply biotin block for 15 minutes at room temperature.

DO NOT RINSE SLIDES WITH BUFFER BEFORE ADDING PRIMARY ANTIBODY.
ONLY WIPE EXCESS BLOCK.

8. Apply primary antibody at a 1:800 dilution. Incubate for 1 hour at room temperature.

Lot # _____ Exp. Date _____

For negative control slides, dilute normal rabbit serum so that its protein concentration matches that of the primary antibody (if necessary). Then make a 1:800 dilution. Apply the negative and incubate for 1 hour at room temperature.

Lot # _____ Date Reconstituted _____

9. Rinse the slides in 2 changes of 1X wash buffer for 5 minutes each.
10. Apply the secondary antibody from Rabbit Elite Kit. Incubate for 30 minutes at room temperature.
11. Rinse the slides in 2 changes of 1X wash buffer for 5 minutes each.
12. Apply the label complex from the Rabbit Elite Kit. Incubate for 30 minutes at room temperature.
13. Rinse the slides in 2 changes of 1X wash buffer for 5 minutes each.
14. Apply the DAB chromogen. Incubate in the dark for 6 minutes at room temperature.
(Add 1 drop of DAB per ml of substrate)
Lot # _____ Exp. Date _____ New Kit: yes / no
15. Rinse the slides in tap water 3 minutes.
16. Counterstain with hematoxylin for 20 seconds.
17. Rinse the slides in tap water until water is clear.
18. Gently agitate slides in 1X wash buffer until the tissues turn blue.
19. Dehydrate through the following solutions:

Solutions	Repetitions	Time
95% Ethanol	1 time	3 minutes
100% Ethanol	3 times	3 minutes
Xylene	2 times	5 minutes

20. Coverslip

Updated 04/13/06