

Detection of COX-1 in Formalin-Fixed, Paraffin Embedded Mouse Tissue

Reagent and Antibody Information

[1X Wash Buffer](#)

[3% Hydrogen Peroxide](#)

[1% BSA Diluent](#)

Distilled Water

[Normal Rabbit IgG – Affinity Purified](#)

[DAB Chromogen](#)

[Hematoxylin](#)

Blocking Serum: Normal Donkey Serum

Jackson ImmunoResearch Laboratories, Inc.

West Grove, PA 19390

www.jacksonimmuno.com

1-800-367-5296

Catalog # 017-000-011

Avidin / Biotin Blocking Kit

Vector Laboratories, Inc.

Burlingame, CA 94010

www.vectorlabs.com

1-800-227-6666

Catalog # SP-2001

Primary Antibody: Rabbit Polyclonal Anti-Mouse COX-1 Antibody

Cayman Chemical

Ann Arbor, MI 48108

www.caymanchem.com

1-800-364-9897

Catalog # 160109

Lot # 0433497-1

Secondary Antibody: Donkey Anti-Rabbit IgG (H+L) Biotin-SP-Conjugated

Jackson ImmunoResearch Laboratories, Inc.

West Grove, PA 19390

www.jacksonimmuno.com

1-800-367-5296

Catalog # 711-065-152

Label Complex: R.T.U. Vectastain Elite ABC Reagent

Vector Laboratories, Inc.

Burlingame, CA 94010

www.vectorlabs.com

1-800-227-6666

Catalog # PK-7100

Staining Procedure

Positive Control Tissue: Male reproductive tract

Stain Localization: Cytoplasmic

1. Deparaffinize and hydrate slides through the following solutions:

Solution	Repetitions	Time
Xylene	2 times	5 minutes
100% Ethanol	2 times	3 minutes
95% Ethanol	2 times	3 minutes
1X Wash Buffer	2 times	5 minutes

2. Quench endogenous peroxidase by placing the slides in 3% hydrogen peroxide for 15 minutes.

3. Rinse the slides in 2 changes of 1X wash buffer for 5 minutes each.

4. Heat-Induced Epitope Retrieval Using The Microwave

Place a full rack of slides into a Tissue Tek® container with 200 ml of distilled water
(Insert blank slides into any empty slots in the rack to ensure even heating of slides)

Microwave for 5 minutes at power level 5.

Cool for 1 minute. (Add more distilled water, if necessary.)

Microwave again for 5 minutes at power level 5. *Temperature Before Cooling Slides* _____

Cool 20 minutes at room temperature.

Rinse the slides in 2 changes of distilled water for 3 minutes each time.

5. Rinse the slides in 2 changes of 1X wash buffer for 5 minutes each.

6. Block with 10% normal donkey serum for 20 minutes at room temperature.

Lot # _____ Date Reconstituted _____

DO NOT RINSE THE SLIDES. CONTINUE TO AVIDIN-BIOTIN BLOCK.

7. Avidin / Biotin Blocking Kit

Lot # _____ Exp. Date _____ New Kit: yes / no

Apply avidin block for 15 minutes at room temperature.

Quick rinse in 1X wash buffer.

Apply biotin block for 15 minutes at room temperature.

DO NOT RINSE SLIDES WITH BUFFER BEFORE ADDING PRIMARY ANTIBODY.
ONLY WIPE EXCESS BLOCK.

8. Apply primary antibody at 1:200 dilutions. Incubate for 1 hour at room temperature.

Lot # _____ Exp. Date _____

For negative control slides, dilute normal rabbit IgG so that it's IgG protein concentration matches that of the primary antibody (if necessary). Then make a 1:200 dilution. If the concentrations can't be matched using this method, the dilution for the negative reagent may need to be adjusted. Apply the

negative and incubate for 1 hour at room temperature.

Lot # _____ Exp. Date _____

9. Rinse the slides in 2 changes of 1X wash buffer for 5 minutes each.

10. Apply the donkey anti-rabbit secondary antibody at a 1:500 dilution. Incubate for 30 minutes at room temperature.

Lot # _____ Date Reconstituted _____

11. Rinse the slides in 2 changes of 1X Wash Buffer for 5 minutes each.

12. Apply the Vectastain R.T.U. Elite Label and incubate for 30 minutes at room temperature.

Exp. Date _____ New Kit: yes / no

13. Rinse the slides in 2 changes of 1X wash buffer for 5 minutes each.

14. Apply the DAB chromogen. Incubate in the dark for 6 minutes at room temperature.

(Add 1 drop of DAB per ml of substrate)

Lot # _____ Exp. Date _____ New Kit: yes / no

15. Rinse the slides in tap water 3 minutes.

16. Counterstain with hematoxylin for 20 seconds.

17. Rinse the slides in tap water until water is clear.

18. Gently agitate slides in 1X wash buffer until the tissues turn blue.

19. Dehydrate through the following solutions:

Solutions	Repetitions	Time
95% Ethanol	1 time	3 minutes
100% Ethanol	3 times	3 minutes
Xylene	2 times	5 minutes

20. Coverslip

Updated 08/15/12